Instruction Manual

Large Capacity Stirrers

EN - English		1
FR - Français		9
ES - Español		18
IT - Italiano		27
DE - Deutsch		36
PT - Português		45
NL - Nederlands		54
NO - Norsk		58
DA - Dansk		62
SV - Svenska		66
FI - Suomi		70
HU - Magyar		74
PL - Polski		78

European Catalogue Numbers:

Model 1000

Euro plug: 444-7115

UK plug: 444-7116

Swiss plug: 444-7117

Model 2000

Euro plug: 444-7118

UK plug: 444-7119

Swiss plug: 444-7120

TABLE OF CONTENTS

Package Contents	1
Warranty	1
Installation	2
Maintenance & Servicing	2
Environmental Conditions	2
Safety Instructions	3
Warning	3
Standards & Regulations	3
Control Panel	4
Specifications	5
Operating Instructions	6-7
Technical Service	8
Troubleshooting	8
EC Declaration of Conformity	82-84

PACKAGE CONTENTS

Model 1000 or Model 2000 Large Capacity Stirrer
 234cm detachable power cord
 Spin bar
 Instruction manual

WARRANTY

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

INSTALLATION

Upon receiving the VWR Large Capacity Stirrer, check to ensure that no damage has occurred in shipment. It is important that any damage that occurred in transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately.

After unpacking, it requires two (2) people to lift the Large Capacity Stirrer, from the bottom of the unit, to place on a level bench, table or floor, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand typical heat produced by the unit. Always place the unit on a sturdy work surface.

The Large Capacity Stirrer is supplied with a power cord that is inserted into the IEC connector on the left side of the unit first, then it can be plugged into a properly grounded outlet. The 120V unit plugs into a 120 volt, 50/60 Hz source. The 230V unit plugs into a 230 volt, 50/60 Hz source.

MAINTENANCE & SERVICING

The Large Capacity Stirrer is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required. It needs no user maintenance beyond keeping the surfaces clean. The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. Spills should be removed promptly. Do not use a cleaning agent or solvent on the front panel or top plate which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR representative.

ENVIRONMENTAL CONDITIONS

Operating Conditions: Indoor use only.

Temperature:	5 to 40°C (41 to 104°F)
Humidity:	20% to 80% relative humidity, non-condensing
Altitude:	0 to 6,562 ft (2000 M) above sea level

Non-Operating Storage:

Temperature:	-20 to 65°C (-4 to 149°F)
Humidity:	20% to 80% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

SAFETY INSTRUCTIONS

Please read the entire instruction manual before operating the Large Capacity Stirrer.

WARNING! DO NOT use the Large Capacity Stirrer in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer, or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift unit by the top plate.

Heavy walled glass or polypropylene container recommended when stirring at high speeds. If spin bar de-couples at higher speeds, thin walled glass containers may break.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the unit or unplug from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly. **DO NOT** immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

STANDARDS & REGULATIONS

VWR international hereby declares under its sole responsibility that the products comply with the requirements of the following directives and associated standards:

Associated EU Directives:

EMC directive	2004/108/EC
LVD directive	2006/95/EC
WEEE directive	2011/19/EU
ROSH directive	2011/65/EU

Safety standards:

IEC 61010-1	Safety requirements for electrical equipment for measurement, control and laboratory use. Part: General Requirements.
IEC 61010-2-051	Part II: Particular requirements for laboratory equipment for mixing and stirring.


UL Std. No. 61010-1

CSA/CAN C22.2 No. 0-M91

CSA/CAN C22.2 No. 61010-1-04

EMC standards:

EN61000-3-2	EN61000-3-3
EN61000-4-2	EN61000-4-3
EN61000-4-4	EN61000-4-5
EN61000-4-6	EN61000-4-11
IEC 61326	

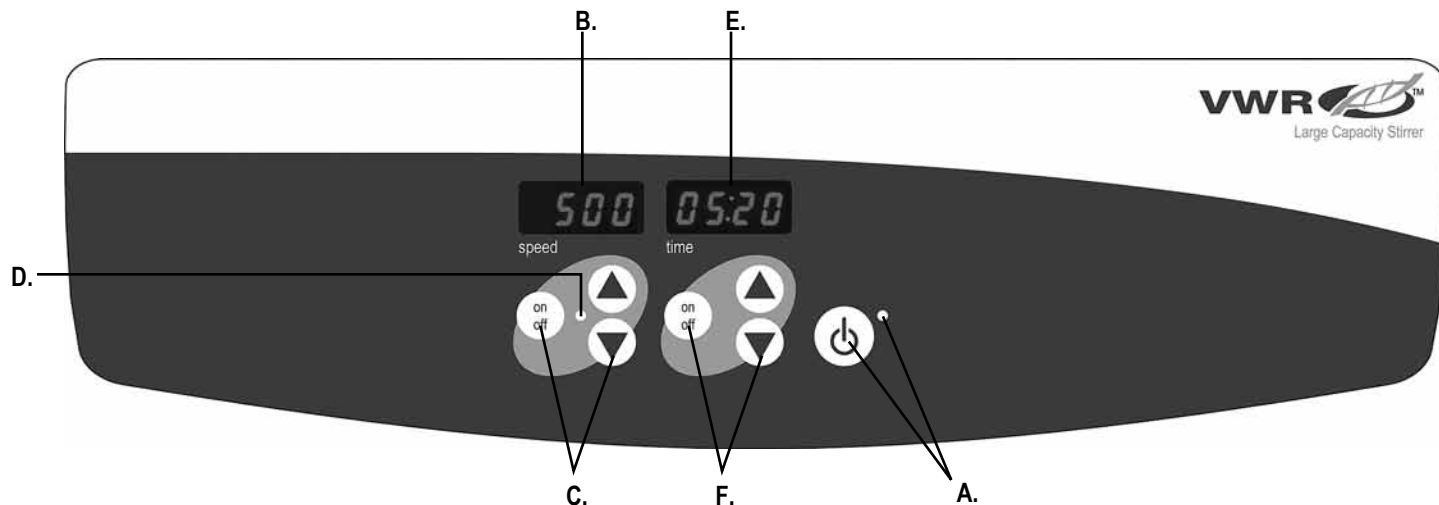CONTROL PANEL

The front panel of the Large Capacity Stirrer contains all the controls and displays needed to operate the unit.

A. Standby button/standby indicator light: The standby indicator light will illuminate when the unit is plugged in. The unit will be in standby mode. Press the standby button to activate the speed and time functions. The standby indicator light will shut off and the speed display and time display will illuminate. Press the standby button again and the unit will once again be in standby mode.

B. Speed display: Displays the speed of the stirrer. **C.** Up/down arrows for set-point control. **D.** The stir indicator light will be illuminated when the unit is stirring.

E. Time display: Displays accumulated time (continuous mode) or how much time is remaining (timed mode). **F.** Up/down arrows for set-point control. On/off button starts/stops the time function. The display range is from 0 to 9,999 minutes in one (1) second increments. The display will indicate minutes and seconds until the timer reaches 99 minutes and 59 seconds (99:59), then the display will automatically display minutes up to 9,999.

SPECIFICATIONS

Overall dimensions (L x W x H):	72,4 x 68 x 13,3 cm
Top plate dimensions (L x W):	54 x 63,5 cm
Electrical (50/60 Hz):	
Model 1000	230 volts 0,5 amps 41 watts
Model 2000	230 volts 0,5 amps 82 watts
Fuses:	5mm x 20mm, 5 amp quick acting
Speed range:	100 to 1800rpm (increased in 10rpm increments)
Speed stability:	±2%
Weight capacity:	193 kg
Volume (H₂O):	
Model 1000	100L
Model 2000	200L
Timer:	digital, 1 second to 9999 minutes (increased in 1 second increments)
Controls:	see page 4
Ship weight:	28,6 kg
European Catalogue Numbers:	
Model 1000	Euro plug: 444-7115 UK plug: 444-7116 Swiss plug: 444-7117
Model 2000	Euro plug: 444-7118 UK plug: 444-7119 Swiss plug: 444-7120

OPERATING INSTRUCTIONS

The Large Capacity Stirrer has been designed for the speed and time functions to work independently of one another. The speed can be re-set without re-setting the timer and the timer can be stopped and started without interrupting the stirring function.

1. Getting ready:

- a. Plug the power cord into a properly ground outlet. The standby indicator light will illuminate, verifying power to the unit.
- b. Press the standby button to move the unit from standby mode. The standby indicator light will turn off and the speed/time displays will illuminate and display the previously used settings.

2. Setting speed:

- a. Press the up/down arrows below the speed display until you reach the desired speed. When you release the arrow button, the unit will blink off and then on indicating the new set speed has been accepted.
- b. Press the on/off button to start the stirring function. The indicator light below the speed display will illuminate to indicate the stirring function is in-use. While accelerating, the indicator light will blink until set-point is reached and then will remain lit until stirring has ceased. The microprocessor controlled ramping feature slowly increases speed until the set-point is reached which helps to avoid splashing and provides excellent low end control. The microprocessor also regulates the stirring speed, sensing your requirements whether stirring an aqueous, viscous, or semi-solid solution.
- c. Speed adjustments can be made without interrupting stirring by using the up/down arrows below the speed display. After the change has been made and you release the arrow button, the unit will blink off and then on indicating the new set speed has been accepted.
- d. To stop the stirring function, press the on/off button below the speed display.

3. Setting time to zero (0:00) and continuous mode: Accumulated time.

- a. and hold the on/off button below of the time display. After three (3) seconds, the display will indicate the previous set time.
- b. Simultaneously press both the up and the down arrows, the display will indicate zero (0:00). The unit time is now set to zero (0:00) minutes. Alternately, you can use the up/down arrows to get to zero (0:00).
- c. Press the on/off button below the time display. The display will indicate accumulated time. The up/down arrows will become inactive. To stop timer, press the on/off button again. **IMPORTANT:** This will **NOT** interrupt the shaking function. Press the on/off button below the speed display to interrupt the shaking function.
- d. To reset, press and hold the on/off button below the time display. After three (3) seconds the display will indicate the previous set time, which was zero (0:00).

4. Setting timed mode: Programmed time.

- a. Press the up/down arrows below the time display until you reach the desired time.
- b. Start this function by pressing the on/off button below the time display. The unit will run for the selected time, the up/down arrows will become inactive while the timer is running. The unit will stop shaking when time display reaches zero (0:00). Four (4) audible beeps will indicate the time down function is complete. The time display will default back to the set time. To repeat this function for the same set time, simply depress the on/off button again.
- c. To interrupt an automatic timing cycle before it is completed, press the on/off button below the time display. The display will flash off and on to indicate the time function is on "hold".

OPERATING INSTRUCTIONS (CONT'D)

IMPORTANT: This will **NOT** interrupt the shaking function. Press the on/off button below the speed display to interrupt the shaking function. Restart the timer by pressing the on/off button below the time display. Unit will continue counting down to zero (0:00). When the display reaches zero (0:00), you will hear the four (4) audible beeps that indicate the count down function is complete and the shaking function will stop.

5. Turning unit off:

- a. To turn the unit off, press the standby button. The speed and time displays will be blank, the standby indicator light will illuminate. The stirrer should be kept in standby mode when not in use. To completely cut off power to the unit, disconnect the power cord from the unit or un-plug from the wall outlet.

OPERATING TIPS

Built-in memory maintains the last used speed and time settings during a power interruption.

The stirrer increases speed at a steady rate until the set-point is reached, if the spin bar is too large or the liquid is too viscous, the stirrer may not reach its set-point or become de-coupled from the units drive magnet. The set-point speed needs to be reduced. The unit is designed to ramp slowly to assure that the spin bar stays coupled.

* The model 2000/200L unit contains a larger spin bar with a center pivoting ring. Ring may need to be replaced after eight (8) hours of use. Contact your VWR representative for replacement spin bars or rings.

TECHNICAL SERVICE

For more information or technical service contact your local VWR representative or visit www.vwr.com.

Web resources: Visit VWR's website at www.vwr.com for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional products information and special offers

TROUBLESHOOTING

Problem	Cause	Solution
E1	Motor locked/failure	Error should not be addressed by the end user. Switch the unit off and contact your VWR Representative for repair.

EN - English		1
FR - Français		9
ES - Español		18
IT - Italiano		27
DE - Deutsch		36
PT - Português		45
NL - Nederlands		54
NO - Norsk		58
DA - Dansk		62
SV - Svenska		66
FI - Suomi		70
HU - Magyar		74
PL - Polski		78

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hereby declare that the products, mentioned in schedule 1 comply with the essential requirements of the following Directives of the European Parliament and council:

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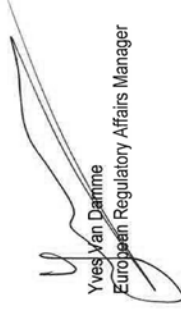
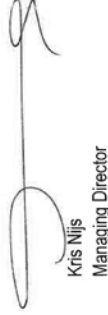
Yves Van Damme
 European Regulatory Affairs Manager

Kris Nijls
 Managing Director

VWR International bvba/sprl Researchpark Haasrode 2020, Geldenaksebaan 464, B-3001 Leuven, Belgium
 Tel.: 016 385 011 • Fax: 016 385 385 • E-mail: info@be.vwr.com

EC Declaration of Conformity

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Schedule 1
Annexe 1
Anhang 1

Article Number	Article description
444-7115	STIRRER LARGE CAPACITY MOD. 100L 230V EU STIRRER LARGE CAPACITY MOD. 100L 230V EU STIRRER LARGE CAPACITY MOD. 100L 230V EU
444-7116	STIRRER LARGE CAPACITY MOD. 100L 230V UK STIRRER LARGE CAPACITY MOD. 100L 230V UK STIRRER LARGE CAPACITY MOD. 100L 230V UK
444-7117	STIRRER LARGE CAPACITY MOD. 100L 230V CH STIRRER LARGE CAPACITY MOD. 100L 230V CH STIRRER LARGE CAPACITY MOD. 100L 230V CH
444-7118	STIRRER LARGE CAPACITY MOD. 200L 230V EU STIRRER LARGE CAPACITY MOD. 200L 230V EU STIRRER LARGE CAPACITY MOD. 200L 230V EU
444-7119	STIRRER LARGE CAPACITY MOD. 200L 230V UK STIRRER LARGE CAPACITY MOD. 200L 230V UK STIRRER LARGE CAPACITY MOD. 200L 230V UK
444-7120	STIRRER LARGE CAPACITY MOD. 200L 230V CH STIRRER LARGE CAPACITY MOD. 200L 230V CH STIRRER LARGE CAPACITY MOD. 200L 230V CH

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Ireland / Northern Ireland

VWR International Ltd /
VWR International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail: sales@ie.vwr.com

Italy

VWR International PBI S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 800 152999/02-40090010
E-mail: info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 200 do 204
Fax: 058 32 38 205
E-mail: info@pl.vwr.com

Portugal

VWR International -
Material de Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43- 3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnologia 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: kundservice@se.vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.ti.
a VWR International Company
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İlmerkezi No.32/1
34896 Pendik - Istanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info@pro-lab.com.tr

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard - Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com

Australia

VWR International, Pty Ltd.
Unit 1/31 Archimedes Place
Murarrie, Queensland 4172
Tel.: 1300 727 696
Fax: 1300 135 123

China

VWR (Shanghai) Co., Ltd
2nd Floor, Building 4,
Lane 998, Halei Rd,
Zhangjiang Hi-tech Park
Shanghai, 201203
Tel.: +86-21-5898 6888
Fax: +86-21-5855 8801
E-mail: info_china@vwr.com

India

VWR Lab Products Private Limited
135/12, Brigade Towers, 2nd Floor
Front wing, Brigade Road,
Bengaluru, India - 560 025
Tel.: +91-80-41117125/26 (Bengaluru)
Tel.: +91-2522-647911/922 (Mumbai)
Fax: +91-80-41117120
E-mail: vwr_india@vwr.com

New Zealand

Global Science - A VWR Company
241 Bush Road
Albany 0632, Auckland
Tel.: 0800 734 100
Fax: 0800 999 002
E-mail: sales@globalscience.co.nz

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

GO TO VWR.COM FOR THE LATEST NEWS, SPECIAL OFFERS AND DETAILS OF YOUR LOCAL VWR DISTRIBUTOR